# REF ASCORBATE HR

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### 1. Positive Interferences (Overestimate Vitamin C/Ascorbate levels)

These compounds consume the ASCOR HR & LR reagent instead of, or in addition to, the ascorbate. The technician will dispense more drops, leading to an overestimation of ascorbate levels.

- **Sulfites:** The most common interference in industry (preservatives, water treatment). They react instantaneously. To avoid this issue, perform steps 3 and 4.
- **Thiosulfates:** Often used to neutralize residual chlorine. They react very strongly with the ASCOR HR & LR reagent.
- **Polyphenols and Tannins:** Found in natural fruit juices (grape, tea, plant extracts). They consume the ASCOR HR & LR reagent more slowly than ascorbate but can skew results by 5 to 10%. If the blue-violet color appears and then disappears, a high concentration of polyphenols may be present.
- **Nitrites:** In an acidic medium (required for the test and managed by the INDIC-ASCORB spoon), they can react and disrupt the reaction equilibrium.

### 2. Negative Interferences (Underestimate Vitamin C/Ascorbate levels)

These compounds destroy the ascorbate before or during the test, or prevent the visualization of the color change.

- **Residual Chlorine (Strong Oxidants):** If the water is chlorinated, the chlorine oxidizes the ascorbate immediately upon its introduction into the water. You will only measure the "surviving" ascorbate.
- **Heavy Metals (Iron III, Copper II):** These ions act as oxidation catalysts. In the presence of oxygen (air), they destroy ascorbate very rapidly.
  - The longer the test takes to perform, the greater the loss.
- **Dissolved Oxygen:** At high pH, atmospheric oxygen oxidizes the ascorbate. This is why the test must be started immediately after sampling.

### 3. Physical and Environmental Interferences

These do not change the chemistry but distort the technician's reading.

- **Temperature:** The powder loses sensitivity above 30-35°C. At high temperatures, the blue-violet color either fails to form or disappears too quickly.
  - **Action:** Always operate on a sample cooled to ambient temperature (20-25°C).
- **pH (Acidity):** If the medium is not acidic enough, the reaction does not proceed correctly. Conversely, if the medium is too acidic, the reaction is also hindered.
  - **Action:** Measure the pH and adjust the amount of INDIC-ASCORB powder if necessary.
- **Natural Water Coloration:**
  - If the sample is very dark (dark fruit juice, wastewater), the transition from clear to blue is invisible.
  - **Solution:** Filter the sample or dilute it with distilled water (and multiply the result by the dilution factor).
- **Intense Light:** The reaction is photosensitive. Titrating in direct sunlight can cause reagent decomposition and skewed results.